



PRECELLYS® EVOLUTION



The most efficient & flexible homogenizer

- Super efficient: DNA, RNA & proteins from any biological sample
- Super strong: powerful 3D motion, up to 10 000rpm
- Super flexible: 6 different volumes of tubes
- Super easy: automatic tube locking system







PRECELLYS® EVOLUTION

Super Homogenizer

The Precellys® Evolution offers optimal robustness, safety, efficiency and flexibility. It is the perfect solution to disrupt samples and get access to DNA, RNA, proteins or small molecules.

Especially designed for L3 laboratories, the Precellys® Evolution is based on 10 years of sample preparation expertise. The instrument's unique figure-8 multi-directional motion provides the same high level of energy and homogenization efficiency to each tube.

The Precellys® Evolution has the capability to process 6 different sizes of tubes, up to 24 tubes simultaneously. Its flexibility and high efficiency gives users the ability to grind, homogenize or lyse any type of soft or hard sample from animal, human, plant or microorganism, in about 30 seconds. Sample are then ready for DNA/RNA analysis, PCR, Electrophoresis, Liquid Chromatography, Western Blot...

Applications







Agronomy



Forensic



Pharmacology




Environment




Microbiology


Precellys® Evolution - Quick User Mode

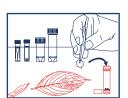

1. Choose the right Precellys Lysing kit (prefilled tubes with beads)

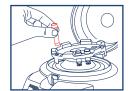

2. Choose the appropriate tube holder and load from 1 to 24 samples

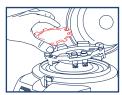

3. Tubes are automatically blocked by the patented tube locking system



4. Select «Use» in the main menu and set up the parameters



5. 3D bead-beating movement for each tube



6. In -30 seconds, your samples are ready for analysis!

🎇 Preserve your sensitive samples & keep your proteins active: ensure the combined yield and integrity of thermo-sensitive molecules by using the Cryolys cooling unit!

Precellys homogenizers range

Discover the full range of products



Precellys® 24



Cryolys® Evolution



Technical specifications

CAPACITY

from 1 to 24 tubes of 2mL/0.5mL from 1 to 12 tubes of 7mL from 1 to 6 tubes of 15mL

from 8 to 96 wells of 300µL from 1 to 6 metal tubes

SPEED RANGE

4 500 rpm to 10 000 rpm (by increment of 100)

PROGRAMS: Unlimited

NUMBER OF CYCLES: up to 10

TIME OF CYCLE: up to 90s

PAUSE BETWEEN CYCLES: 120s

L: 378mm, W: 515mm, H: 400mm (632mm with opened lid)

WEIGHT: 27,2 kg

NORMS: CE, FCC part 15

ELECTRICITY

90 - 250VAC / 50-60 Hz power:1 kVA

NOISE < 70dB

CATALOG NUMBER P000062-PEVO0-A